

BioLane HTI 16V



BioLane HTI 16V

Automated In Situ Hybridization and Immunohistochemistry



- History
- Applications
- Hardware
- Accessories
- Software
- BioLane HTI 16V for Western blot processing




- Developed in 2001 by H&H in cooperation with JOBO as an automated system for (Western) blot staining
- *In situ* application officially launched in 2003 with 48-well plates; especially popular in zebra fish community
- Introduction of individual baskets 2004
- Development of the slide tub 2005
- June 2008, Intavis and H&H instrument joined together
- In 2008, development of the new BioLane HTI 16V, an advanced version of the old BioLane HTI
- Totally installed units about 170, exclusively Western 60-70 (of these for prionics test ~35)

BioLane HTI 16V – Applications


BioLane HTI 16V



In situ hybridization

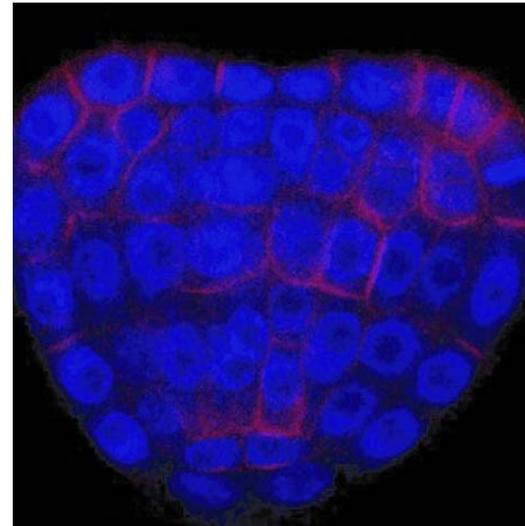

Immunohistochemistry

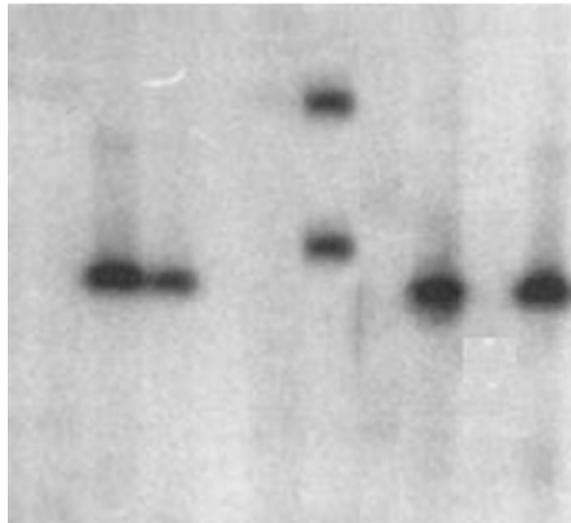

Western blot processing

Hardware – Two systems in one

BioLane HTI 16V



Hardware – Two systems in one


BioLane HTI 16V



- Two independent incubation units
 - heating/cooling
 - shaking

Hardware – Two systems in one


BioLane HTI 16V

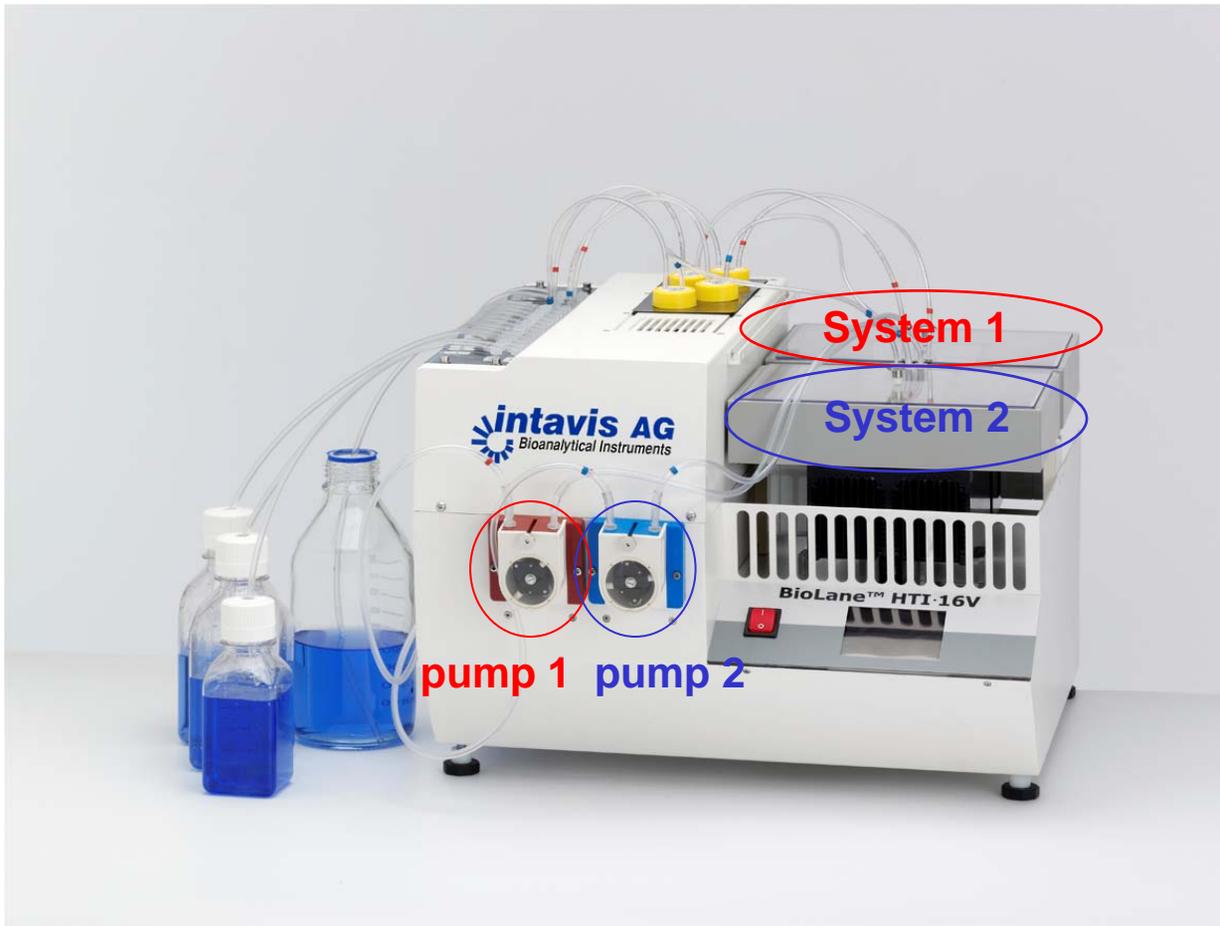

- Two independent incubation units
 - heating/cooling
 - shaking
- Two independent peristaltic pumps
 - for buffer delivery and recovery

Hardware – Two systems in one

BioLane HTI 16V

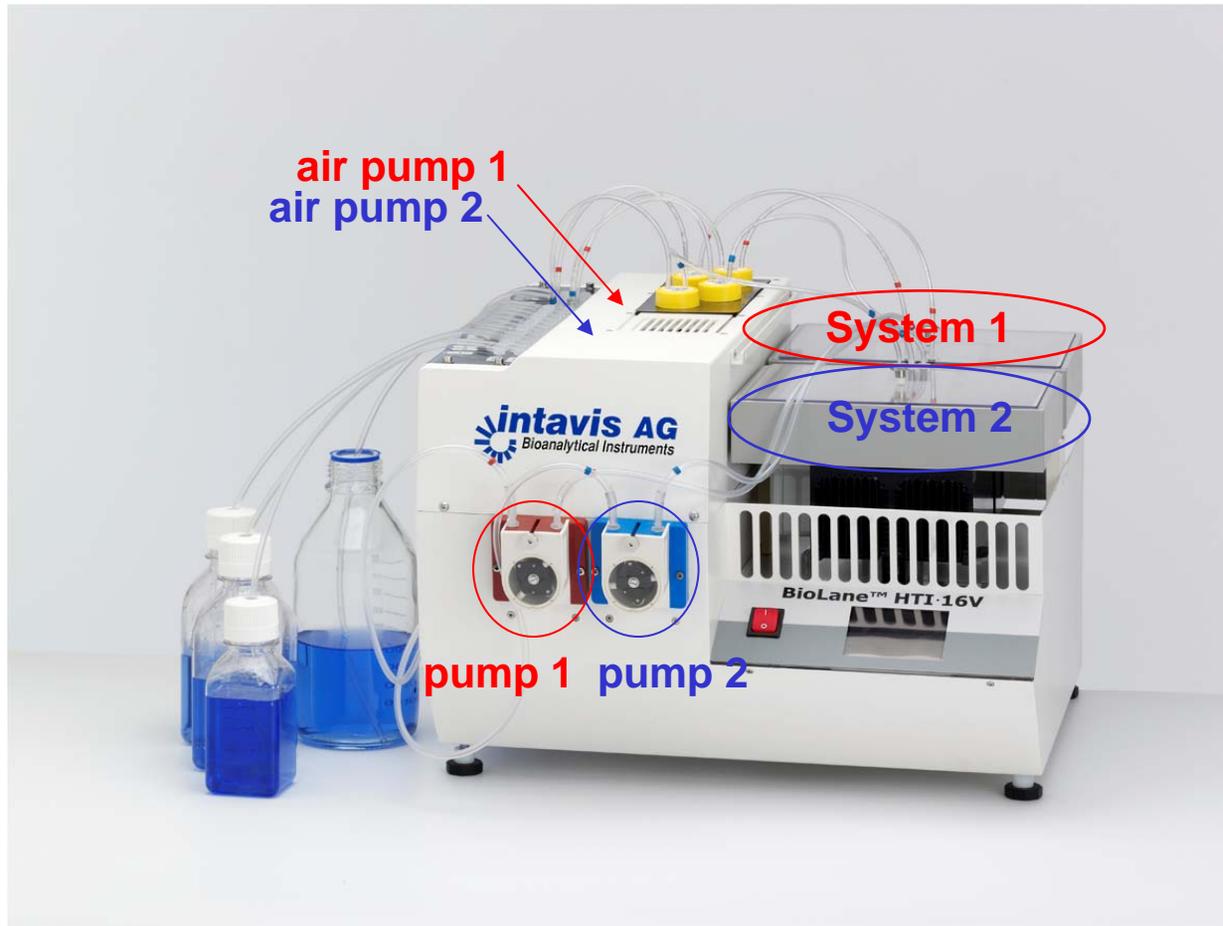

- Two independent incubation units
 - heating/cooling
 - shaking
- Two independent peristaltic pumps
 - for buffer delivery and recovery
- **Two independent air pumps**
 - for antibody delivery and recovery

Hardware – Two systems in one

BioLane HTI 16V



- Two independent incubation units
 - heating/cooling
 - shaking
- Two independent peristaltic pumps
 - for buffer delivery and recovery
- Two independent air pumps
 - for antibody delivery and recovery
 - **Cooled positions**
 - 2x2 positions for 50 ml Falcons or 125ml vials
 - individual tube paths

Hardware – 16 way valve

BioLane HTI 16V



16 way valve

- One 16 way valve for buffer selection
 - 2x 14 positions for buffers
 - 2x 2 positions for waste
 - independent connections for both systems




- Operated by touch screen PDA
 - running the instrument easy by PDA
 - LAN connection for online monitoring
 - programming via external PC or PC in the network via remote control



- Tray plus accessories for application
 - + Medium baskets for ISH/IHC up to 40 samples/unit
 - + Large baskets for ISH/IHC of larger specimens up to 24 samples/unit
 - + plates for ISH/IHC screen up to 48 samples/unit
 - + Slides for washing ISH/IHC of sections up to 25/unit
 - Only trays for Western Blot up to 4 blots/unit, depending on size
 - + 2D incubation unit and trays for 2D-Western, up to two/unit

Accessories – trays

BioLane HTI 16V

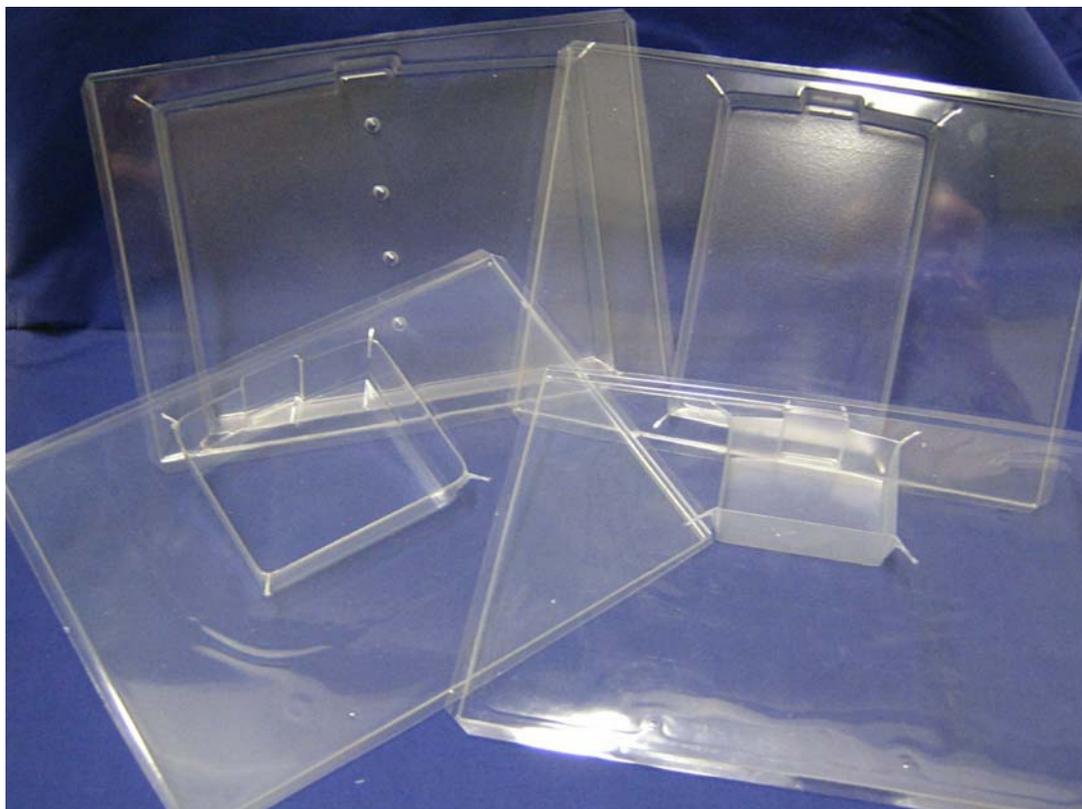

- **4 different tray sizes**

- Extra small 12-16ml
- Small 25-30ml
- Medium 50-60ml
- Double medium 100-120 ml

Accessories – incubation baskets

BioLane HTI 16V



Extra small:

- Up to 8 medium

Small:

- Up to 20 medium

Medium:

- Up to 40 medium



Extra small:

- Up to 6 large

Small:


- Up to 12 large

Medium:

- Up to 24 large

Accessories – medium baskets hybridization


BioLane HTI 16V

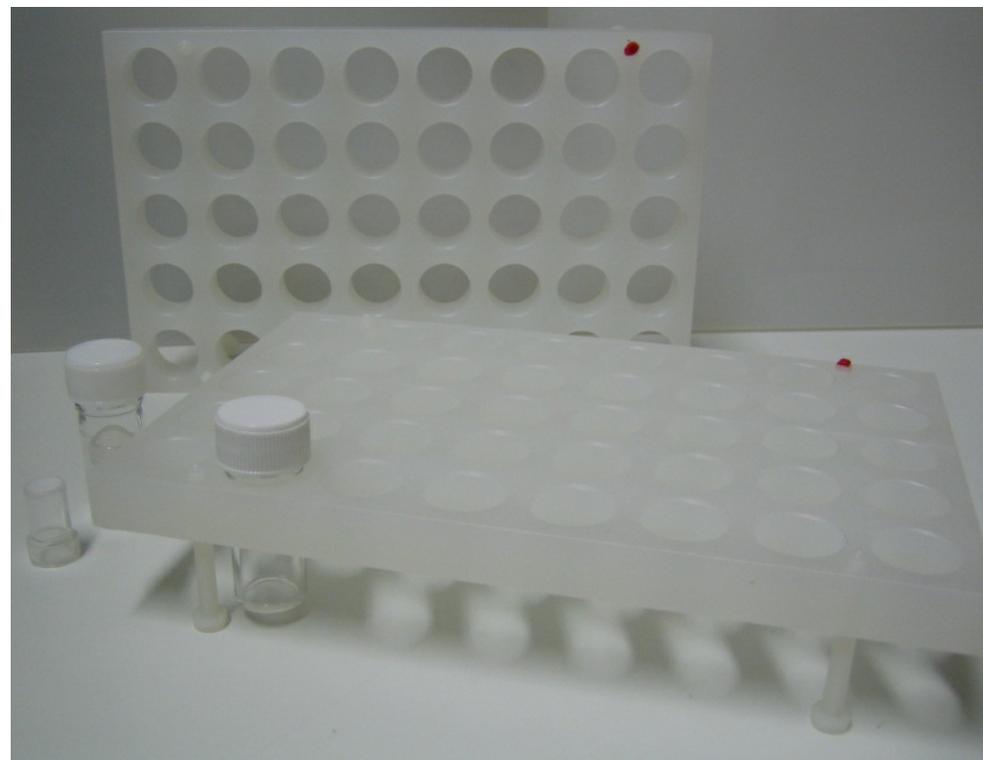

Hybridization in vials

Accessories – large baskets hybridization


BioLane HTI 16V

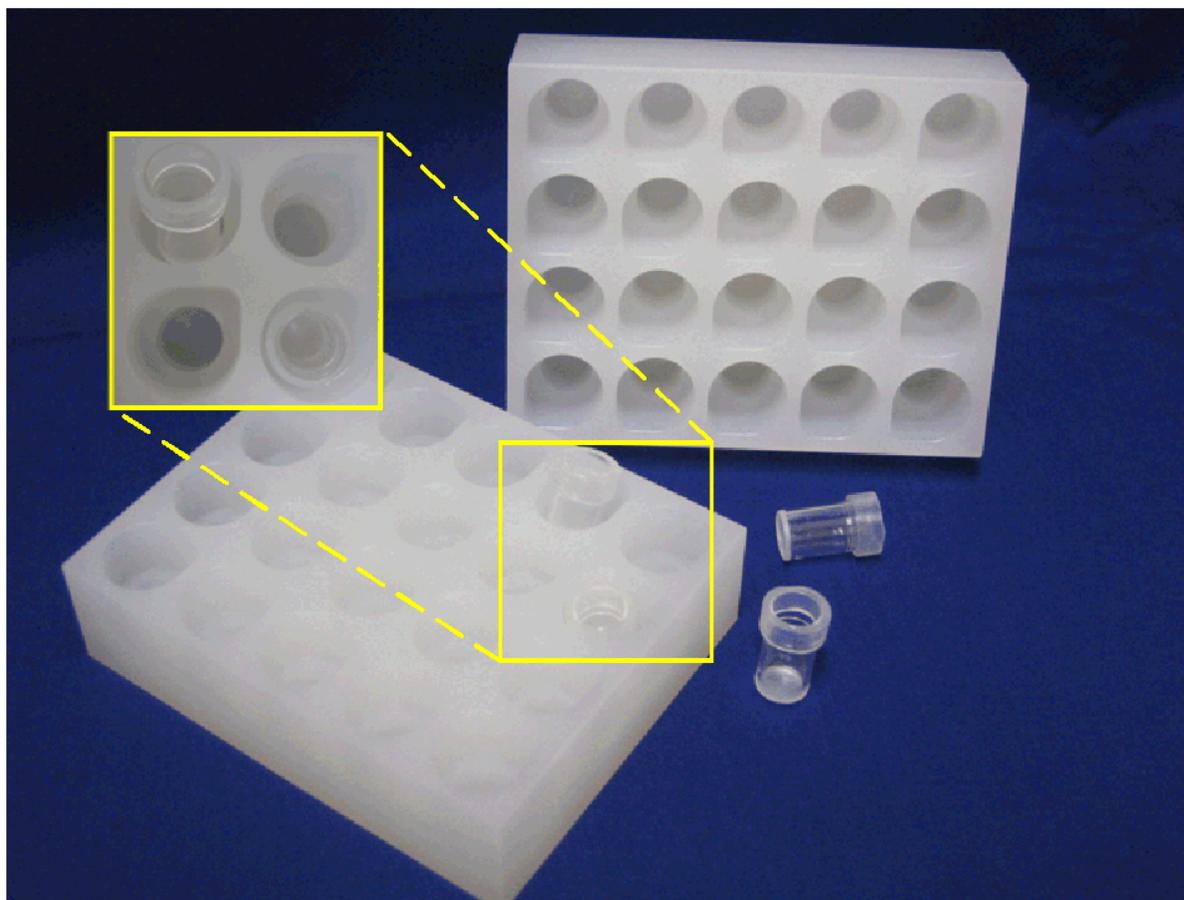

Hybridization in plate

Accessories – plate and plate incubation



BioLane HTI 16V

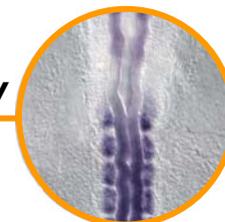


Only for medium tray: 48 medium in plate



Accessories – slide incubation

BioLane HTI 16V

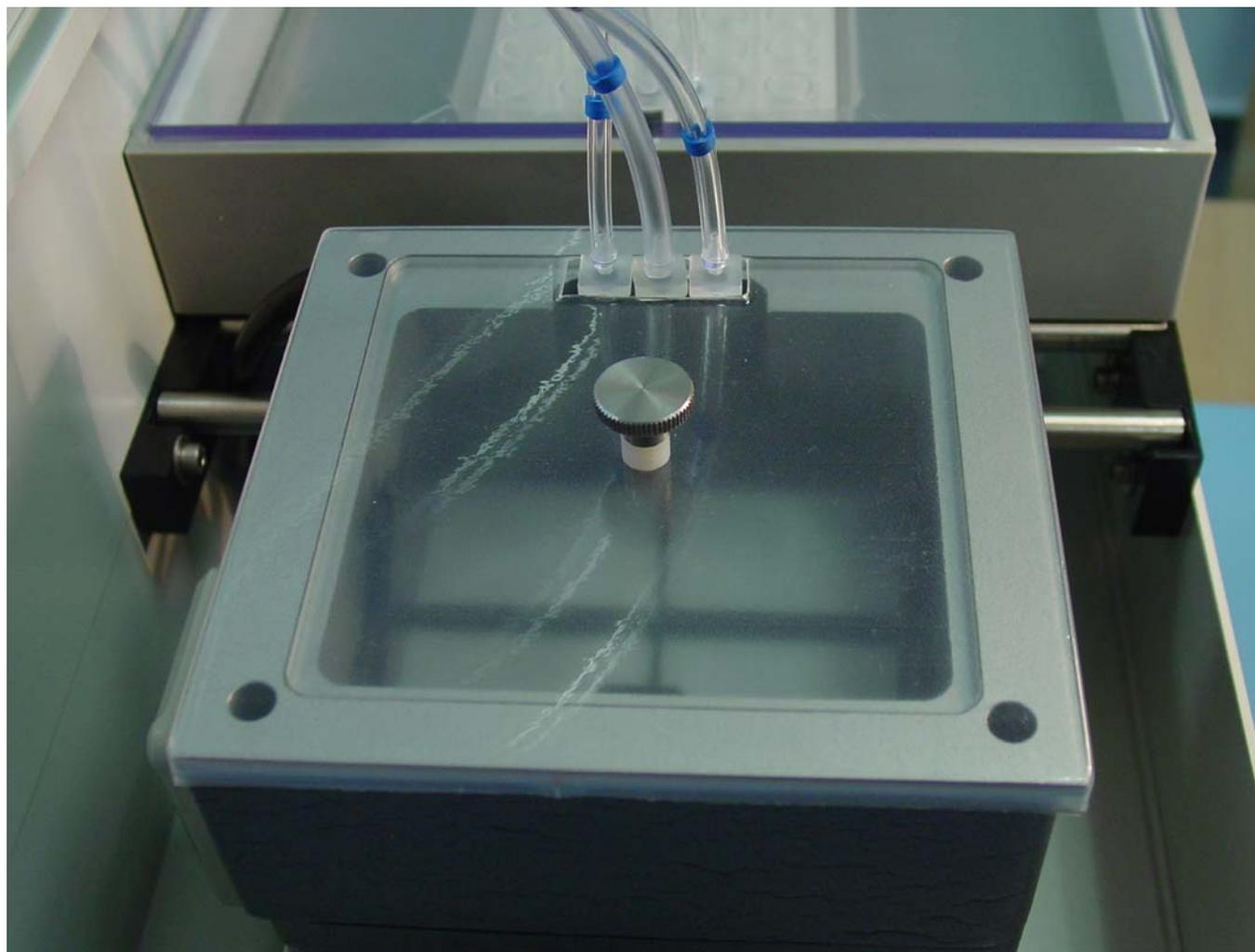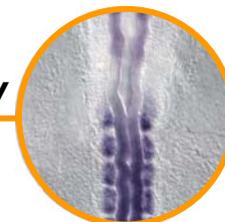

- Only washing
- Up to 25 slides/ incubation unit

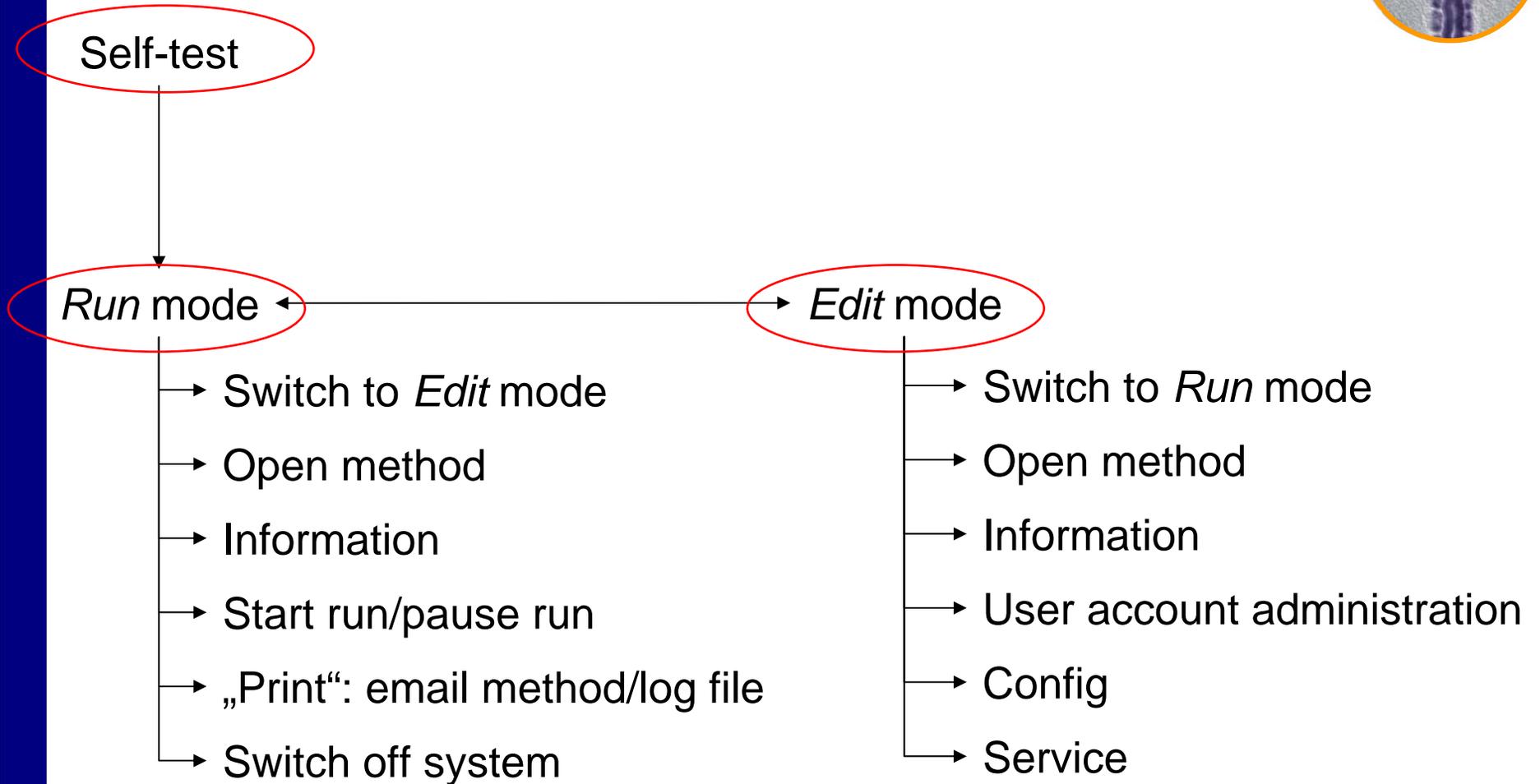

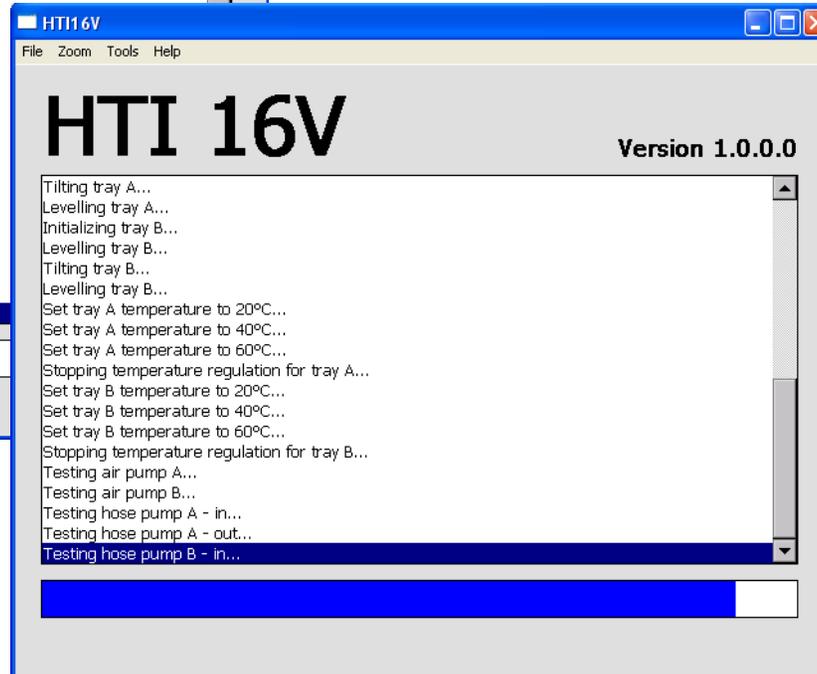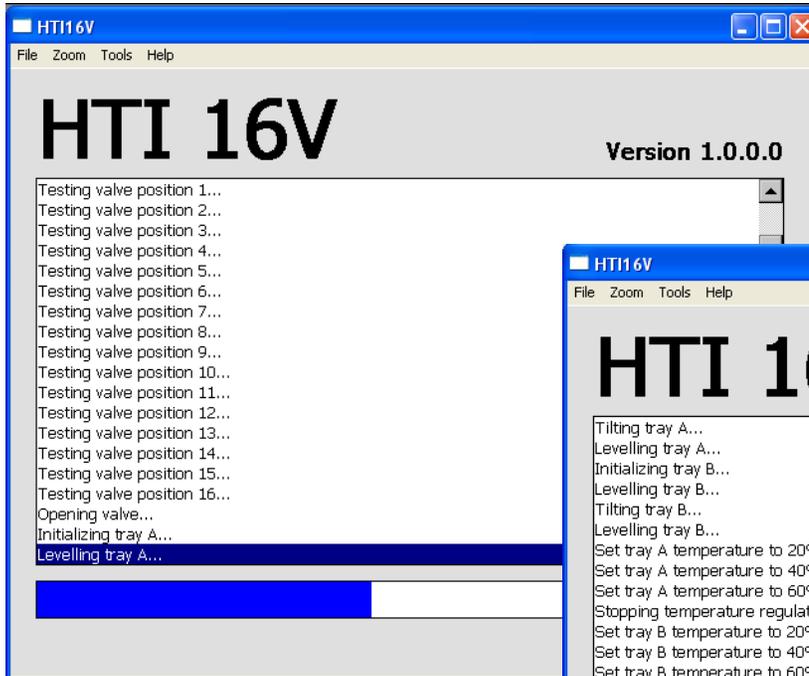

Software – The Run mode

BioLane HTI 16V



HTI16V

File Zoom Tools Help

1:50:06 PM Antibody: 22.0 °C Heater: n/a Current User: -

1/23/2009

Owner: Lab
Name: In Situ
Current Tray: Small

System 1
Idle
23.1 °C
1.23:09:30

System 2
Idle
23.5 °C
1.23:09:30

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	...
Heat			Sol I			Sol I									
Agitation [1/min]: 0			Agitation [1/min]: 25			Agitation [1/min]: 25									
Temp. [°C]: 60.0			Temp. [°C]: 60.0			Temp. [°C]: 60.0									
Pause after: No			Pause after: No			Pause after: No									
Pause before: No			Pause before: No			Pause before: No									
Time [h:m:s]: 00:05:00			Time [h:m:s]: 00:15:00			Time [h:m:s]: 00:15:00									
Drain: Valve B			Drain: Valve B			Drain: Valve B									
00:00:00			00:00:00			00:00:00									

1. Status line
2. TCP/IP button to connect to a network
3. System information
4. Method navigation arrows
5. Method step numbers
6. Red system (system 1)
7. Blue system (system 2)
8. Combi step
9. Icon bar

Software –The Edit mode

BioLane HTI 16V



HTI16V

File Zoom Tools Help

Steps: [Icons] Current User: Administrator

Owner: Lab Base Tray: Extra Small

Name: In Situ Current Tray: Small


1 2 3 4 5 7 8 9 10 11 12 13 14 15 16 17 18 19 20 ...

Heat	Sol I	Sol I	Sol I
Source: Valve 1	Source: Valve 1	Source: Valve 1	Source: Valve 1
Vol.[ml]: 16.0 (Base)	Vol.[ml]: 16.0 (Base)	Vol.[ml]: 16.0 (Base)	Vol.[ml]: 16.0 (Base)
Agitation [1/min]: 0	Agitation [1/min]: 25	Agitation [1/min]: 25	Agitation [1/min]: 25
Temp. [°C]: 60.0	Temp. [°C]: 60.0	Temp. [°C]: 60.0	Temp. [°C]: 60.0
Pause after: No	Pause after: No	Pause after: No	Pause after: No
Pause before: No	Pause before: No	Pause before: No	Pause before: No
Time [h:m:s]: 00:05:00	Time [h:m:s]: 00:15:00	Time [h:m:s]: 00:15:00	Time [h:m:s]: 00:15:00
Drain: Valve B	Drain: Valve B	Drain: Valve B	Drain: Valve B
00:00:00	00:00:00	00:00:00	00:00:00

00:00:00 00:00:00 00:00:00 00:00:00

[Play] [Folder] [Info] [People] [Gear] [Wrench]

1. Status line
2. Small icon bar to edit programs
3. Program information
4. Method navigation arrows
5. Method step numbers
6. Inpump step
7. Incubation step
8. Outpump step
9. Icon bar

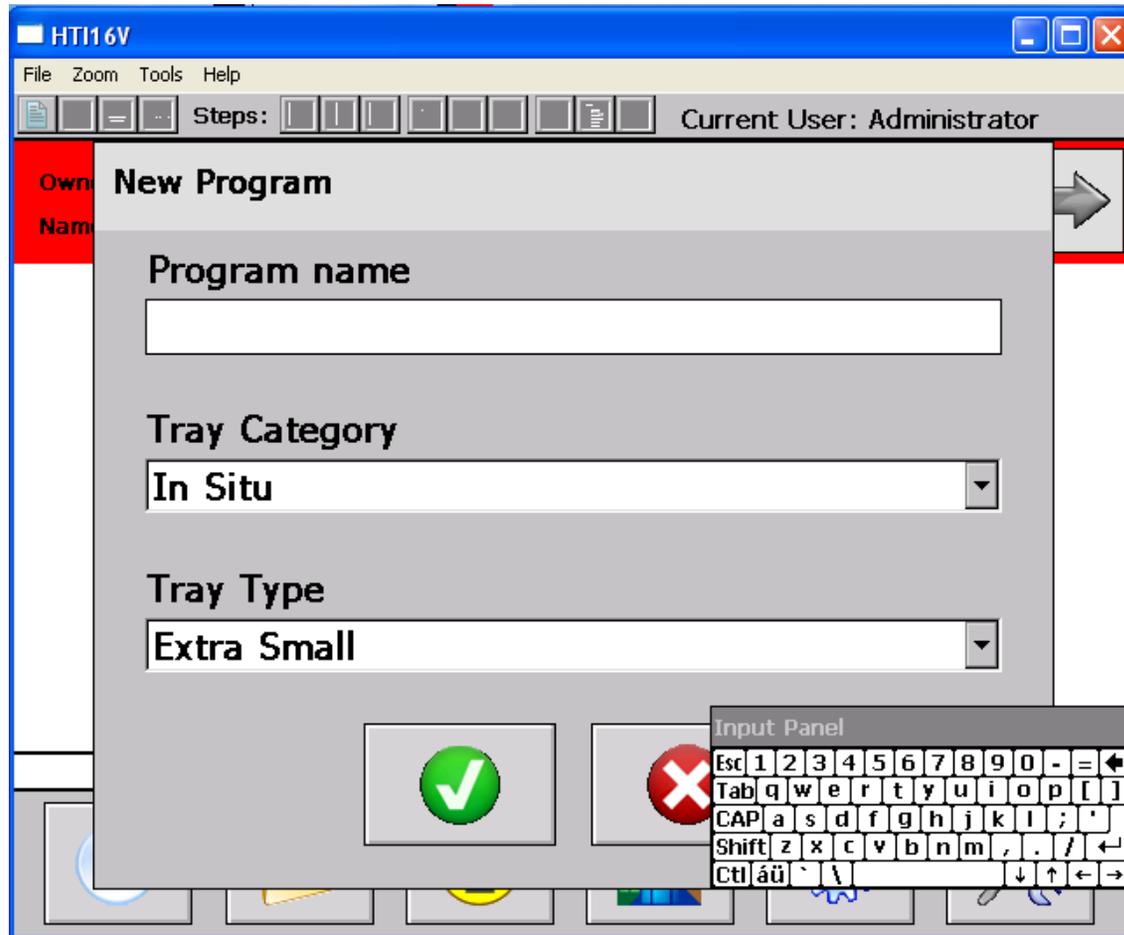

- Program name
- Tray category
- Base tray type

Software – Steps

BioLane HTI 16V



HTI16V

File Zoom Tools Help

Steps: [Navigation icons] Current User: Administrator

Owner: Administrator Base Tray: Extra Small
Name: In Situ Day 1 Current Tray: Extra Small

1 2 [3] 4

PBST	PBST	Blocking	PBST
Source: Valve 1 Vol.[ml]: 20.0 (Base)	Source: Valve 1 Vol.[ml]: 20.0 (Base)	Source: Valve 1 Vol.[ml]: 20.0 (Base)	Source: Valve 1 Vol.[ml]: 20.0 (Base)
Agitation [1/min]: 25 Temp. [°C]: 22.0 Pause after: No Pause before: No Time [h:m:s]: 00:05:00	Agitation [1/min]: 25 Temp. [°C]: 22.0 Pause after: No Pause before: No Time [h:m:s]: 00:05:00	Agitation [1/min]: 25 Temp. [°C]: 4.0 Pause after: No Pause before: No Time [h:m:s]: 00:05:00	Agitation [1/min]: 25 Temp. [°C]: 22.0 Pause after: No Pause before: No Time [h:m:s]: 00:05:00
Drain: Valve B 00:00:00	Drain: Valve B 00:00:00	Drain: Valve B 00:00:00	Drain: Valve B 00:00:00

[Control icons]

- Step name
- Inpump step
- Incubation step **Combistep**
- Outpump step
- Progress bar



HTI16V

File Zoom Tools Help

Information

Volume Time

Position	Volume
Valve 1	142.8
Valve 2	35.7
Valve 3	142.8
Valve 4	35.7
Valve 5	178.4
Valve 6	856.6
Valve 7	107.1
Valve 8	0.0
Valve 9	0.0
Valve 10	0.0
Valve 11	0.0

Tray Category: In Situ
 Base Tray Type: Extra Small
 Current Tray Type: Small

✓ ✗

Volume information

HTI16V

File Zoom Tools Help

Information

Volume Time

Total Time: 47:09:30

Step #	Duration [hh:mm:ss]	1 to...	2 to...	3 to...	4 to...	5 to...	6 to...
1	00:05:30						
2	00:16:00	00:05:30					
3	00:16:00	00:21:30	00:16:00				
4	00:16:00	00:37:30	00:32:00	00:16:00			
5	00:16:00	00:53:30	00:48:00	00:32:00	00:16:00		
6	00:11:00	01:09:30	01:04:00	00:48:00	00:32:00	00:16:00	
7	00:06:00	01:20:30	01:15:00	00:59:00	00:43:00	00:27:00	00:
8	00:06:00	01:26:30	01:21:00	01:05:00	00:49:00	00:33:00	00:
9	00:06:00	01:32:30	01:27:00	01:11:00	00:55:00	00:39:00	00:

✓ ✗

Runtime information


Software – Online monitoring and remote control BioLane HTI 16V




- System can be directly connected to a PC by a crossed cable or to a LAN network using TCP/IP
- Via remote control, you can directly operate the system or monitor progress online
- Reminders can be send by email when program finished
- Protocols can be sent by email

BioLane HTI 16V – Applications


BioLane HTI 16V

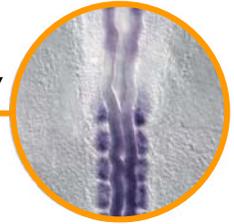

In situ hybridization

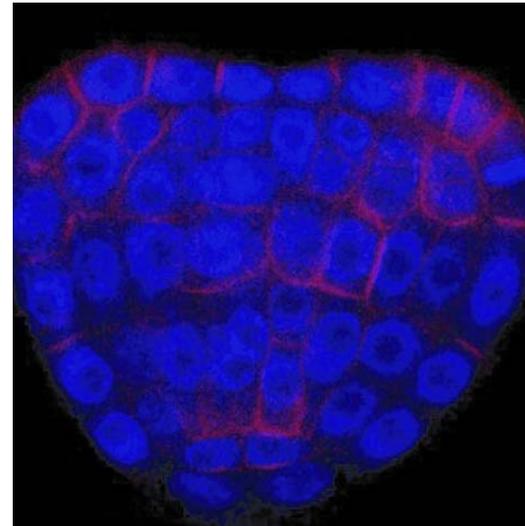

Immunohistochemistry

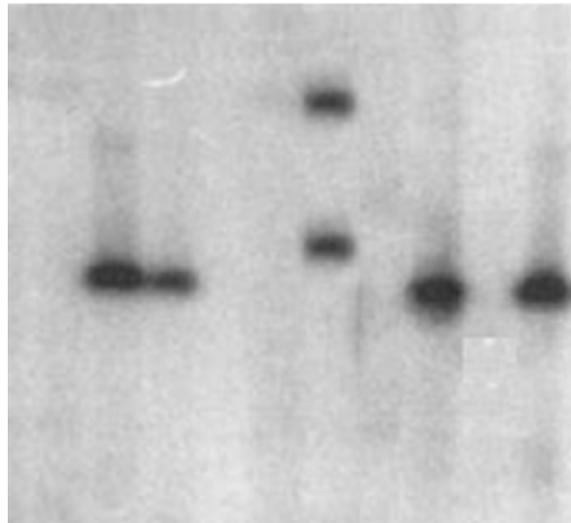

Western blot processing



- Separation of a protein mixture by SDS-gel electrophoresis
- Transfer to and immobilization on membrane
- Processing for detection

typically

3x washing
primary antibody
3x washing
secondary antibody
3x washing
Detection

Western Blotting

Western Blot
Processing

Can be automated by
the BioLane HTI 16V

Western Blot Processing – Why BioLane HTI 16V

BioLane HTI 16V



- Improves results
 - ◆ Independent tube paths for buffers, primary and secondary AB
 - ◆ Antibody binding easy to optimize due to temperature control
 - ◆ Luffing, no round shaking for max liquid exchange
 - ◆ Fluorescent antibodies can be used due to non-transparent lid
- Increases reproducibility:
 - ◆ Time, volume, temperature reproducibly the same, independent from environmental conditions
 - ◆ Standardization even for untrained personal
 - ◆ Automated GLP reports possible
 - ◆ Better comparability due to standardized protocols. Better initial data → better quantification results (especially 2D gels)
- Saves time
 - ◆ System can work o/n as well
 - ◆ Higher temperature allows shortening of incubation times
- Can handle and recover antibodies
 - ◆ Uninterrupted cold chain
 - ◆ Vial does not have to be changed
 - ◆ Higher consumption, but can be recovered and reused

Western Blot Processing – Competition

BioLane HTI 16V



Competitors

- LI-COR Blot Washer
- Stovall Blot Washer
- Freedom Rocker
and others...

→ Simple tumblers with pumping system

Compared to BioLane HTI 16V

- No significant improvement of results
 - ◆ Usually no independent tube path for antibody
 - ◆ No easy optimization of binding conditions
- Limited increase in reproducibility
 - ◆ Usually no heating/cooling of incubation units
- Time saving limited
 - ◆ Usually no automated application of antibody
- Unsufficient antibody handling
 - ◆ No uninterrupted cold chain or recovery



Li-COR Blot washer


BUT: cheap!

Western Blot Processing – Potential customers

BioLane HTI 16V



- Prionics (BSE testing)
- Clinical applications, e.g. high safety labs (S3+/L3+)
- 2D-Western
- Industrial (GLP-reports)

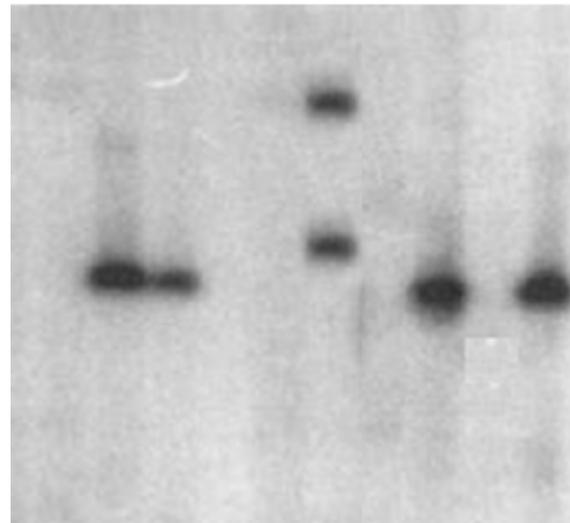

Western blot